Infrico medcare

We care
about your
success

PHARMACARE

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INTERIOR FORCED VENTILATION ENSURES MINIMUM TEMPERATURE VARIATIONS AND ADEQUATE HUMIDITY FOR THE PRESERVATION OF MEDICINES AND THEIR PACKAGING.

PRODUCT STABILITY RANGE OF APPROXIMATELY $\pm 1^{\circ}\text{C}$ AND PRODUCT UNIFORMITY TEMPERATURE LESS THAN $\pm 2^{\circ}\text{C}$.

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FORCED-AIR EVAPORATOR

- Uniform cooling.
- Quick recovery after door opening.
- Better internal humidity control.

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We have our own laboratory with a climatic chamber, dedicated exclusively to test our equipment.

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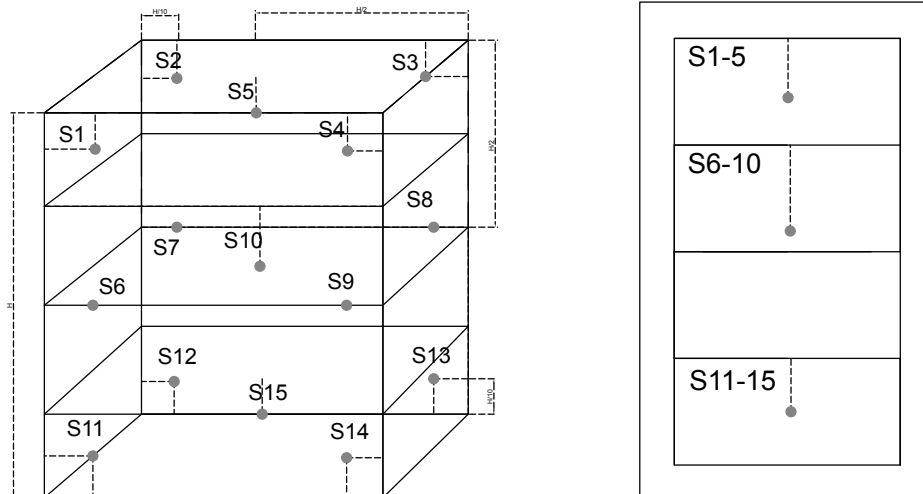
INFRICO MEDCARE TESTS ITS EQUIPMENT IN DIFFERENT CLIMATE CONDITIONS AND PERFORMS A MAPPING OF TEMPERATURES IN THE AIR AND OF THE PRODUCT THROUGHOUT THE INTERIOR OF THE CHAMBER. ALL THIS, IS TO GUARANTEE THE HOMOGENEITY OF THE TEMPERATURE WITHIN THE ESTABLISHED RANGES.


The devices used to make the tests are 9 probes situated according DIN 13277.

- Probe in each shelf—control of the air temperature in the different levels of the equipment
- Probe immersed in liquid—control of the product temperature -
- Probe at the air outlet—control of the temperature of the air outlet, at different levels
- Probe in exchanging air—control of the air temperature in the entrance to the evaporator

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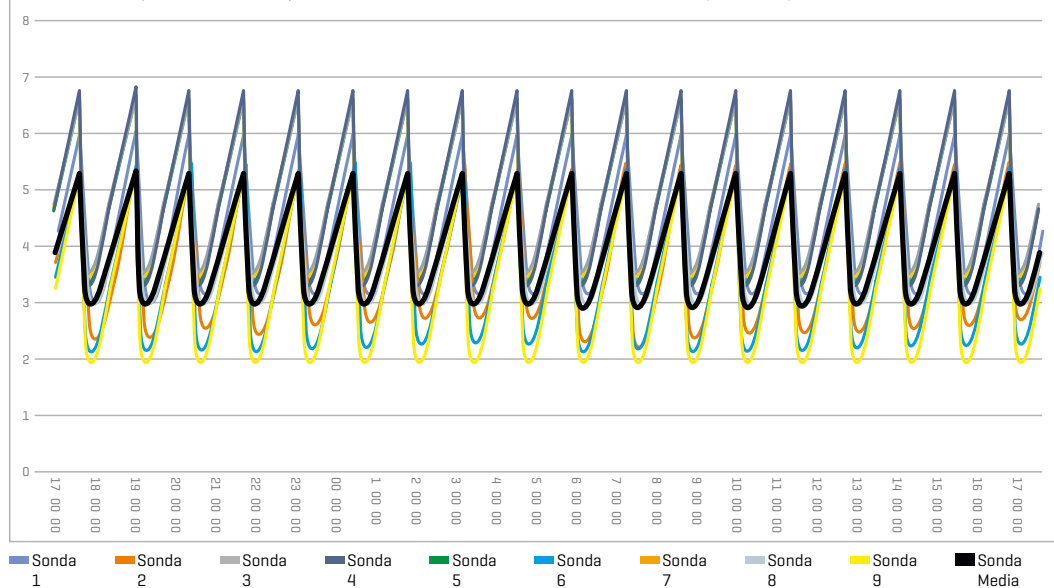EASY AND INTUITIVE - touch screen
 PRECISE CONTROL - NTC probes
 STABILITY AND UNIFORMITY - product stability range of approximately $\pm 1^{\circ}\text{C}$ and product uniformity temperature less than $\pm 2^{\circ}\text{C}$
 GREAT SIZE DISPLAY - good visibility and 0.1°C accuracy
 3 TYPES OF OPTICAL AND ACOUSTIC ALARMS - temperature rise or fall, probe failure and open door
 REMOTE ALARM SIGNAL - Potential free contact
 MULTIPLE OPTIONAL FUNCTIONS - USB, Datalogger Bluetooth or Wifi, Battery, etc.

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USB
 Easy download of data logs and alarms.

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CONNECTIVITY
 Possibility of monitoring via Bluetooth or WIFI.

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BATTERY
 Allows the maintenance of temperature and alarms register in case of a power failure.

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DIFFERENT LANGUAGES SELECTION OPTION
 Spanish, English, Italian, French, German, Turkish and Polish.

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RS485 COMMUNICATION PORT MODBUS
 Connects the system with PC or network.

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REMOTE ALARM SIGNAL
 Potential free contact, allows the transfer of warning signals to an external system.

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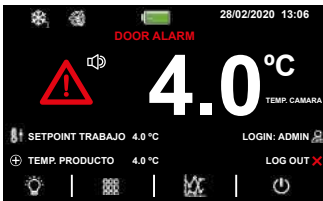
DISPLAY:
Accuracy of 0.1°C.

PRECISE CONTROL:
PT1000 probes

STABILITY AND UNIFORMITY:
Product stability range of approximately $\pm 1^\circ\text{C}$ and product uniformity temperature less than $\pm 2^\circ$.

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
VARIETY OF OPTICAL AND ACOUSTIC ALARMS
Temperature rise or fall, power failure, low battery, open door, compressor maintenance, dirty condenser...


DATALOGGER
Registro de temperaturas e histórico de alarmas

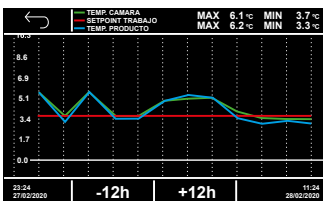DATALOGGER
Temperature log and alarm history

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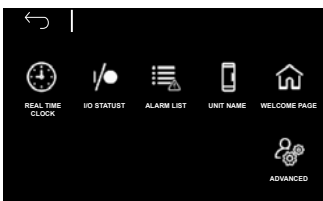
DETAILED EQUIPMENT INFORMATION
5 inches Graphic interface indicating maximum and minimum temperatures

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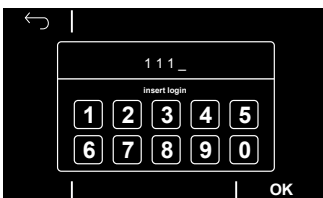
CONFIGURATION ASSISTANT
Welcome page for quick and easy configuration by choosing the main controller parameters

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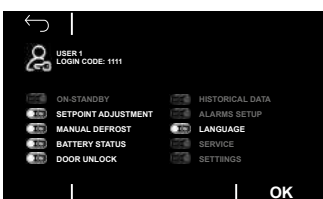LOCK CODE
Lock code to access the control system

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CONFIGURABLE USERS
Permission management to different users

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The sector of pharmacies, hospitals and medical practices is especially demanding in terms of temperature stability. Infrico Medcare in its Pharma CARE and Pharma CARE PLUS product lines has incorporated a series of specific characteristics that guarantee the safety in the conservation of medicines and vaccines 24 hours a day and 365 days a year. This is crucial for our customers.

All our refrigerators of the Pharma CARE PLUS and Pharma CARE PLUS lines are designed in compliance with DIN 58345.

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According to DIN 13277 the devices must comply with the following requirements:

Operating temperature range from +2°C to +8°C.

Interior room temperature of the installation site is accepted between +10°C and +35°C.

Interior is easy to clean and disinfect; water resistant and with corrosion protection.

Mechanical capacity load of uniformly mounted parts of 100 kg/m².

Doors fitted with locks.

Acoustic emission below 60 dB [A]

Power failure alarm for a minimum of 12 h.

Optical-acoustic alarm in case of temperature deviation.

Potential free alarm contact.

Safety device to prevent the temperature going below +2°C

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Using INVERTER technology in our equipment, we achieve greater temperature stability, lower noise level and outstanding energy efficiency.

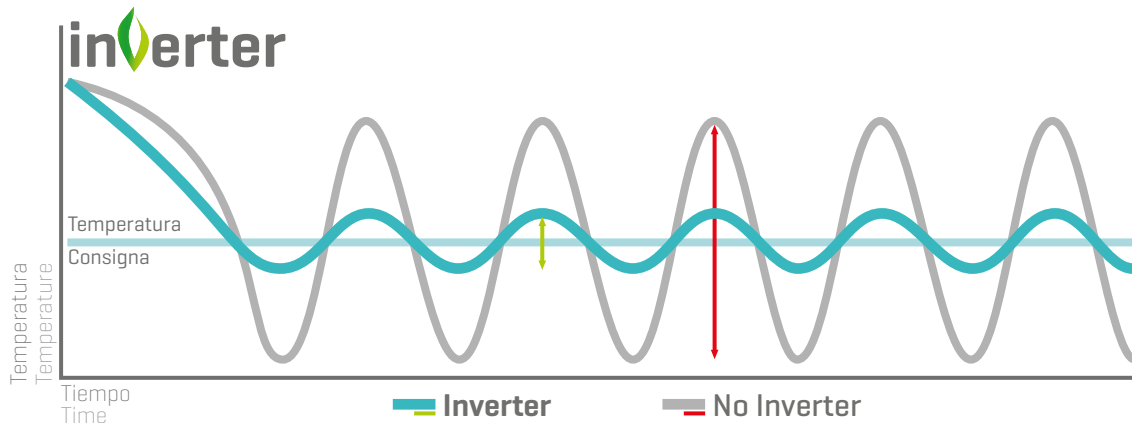
DESTACADA EFICIENCIA ENERGÉTICA
OUTSTANDING ENERGY EFFICIENCY

MENOR CONSUMO ENERGÉTICO
LOWER ENERGY CONSUMPTION

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

Conventional non-INVERTER equipment, the compressor operates at maximum speed until reaching the setpoint temperature, and once it is reached, it stops, repeating successively with continuous starts and stops, causing greater energy consumption and lower temperature stability.

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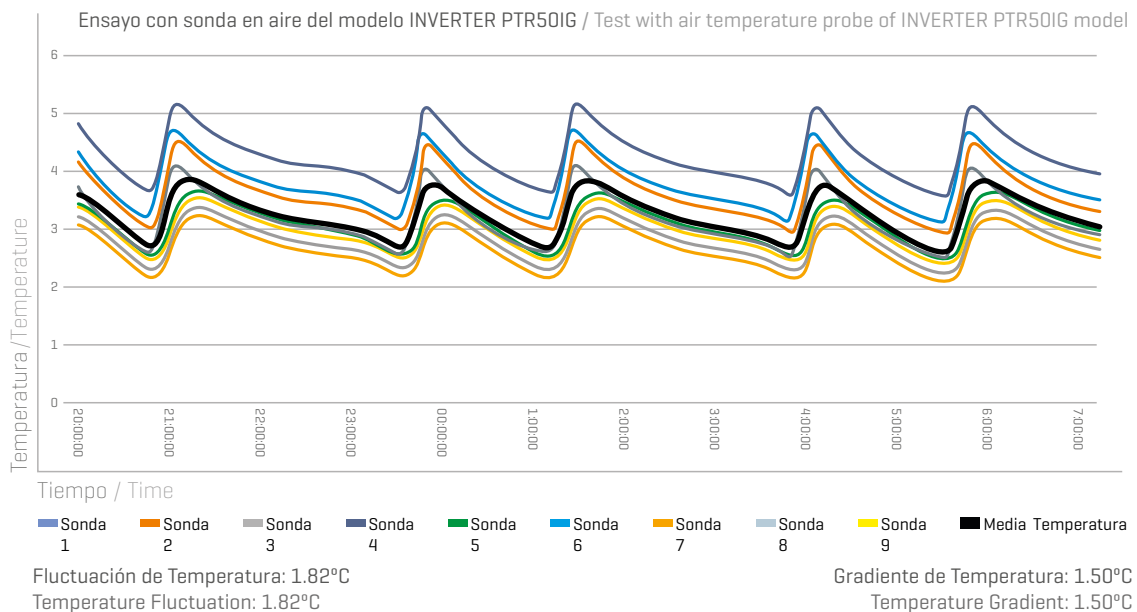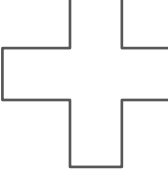In equipment with INVERTER technology, once the compressor starts operating, it adjusts the operational speed to the setpoint temperature, avoiding the starts and stops of the compressor that increase energy consumption and make greater the oscillation of the equipment temperature.

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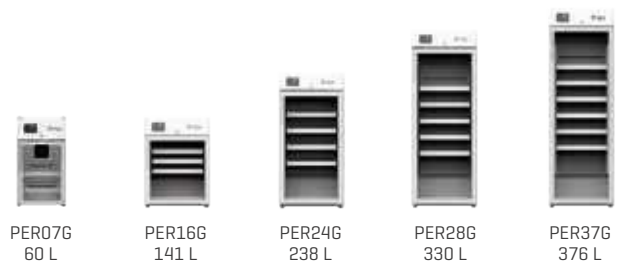
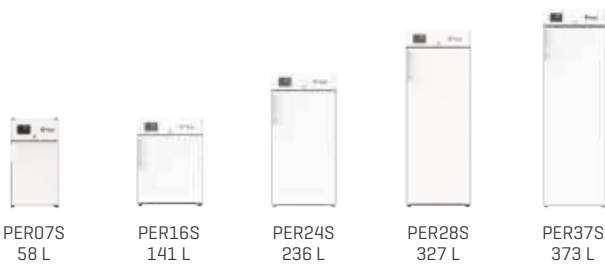
OUR EXTENSIVE EXPERIENCE AS A WORLDWIDE MANUFACTURER OF COOLING EQUIPMENT AND THE SPECIFICATIONS OR QUALITY REQUIREMENTS OF DIFFERENT MARKETS THAT WE COMPLY WITH, ENSURES YOU CAN RELY ON INFRICO MEDCARE.

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PHARMA CARE SERIES

Pharma CARE SERIES offers pharmaceutical refrigerators in different sizes and storage capacities, in order to fit all possible space requirements and needs. Models range from countertop, underbench or upright versions which are available with model variety from 58 to 376 litres. These models are designed according to **DIN 58435** standard. Solid or glass doors are available in the whole range.

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
PHARMA CARE PLUS SERIES

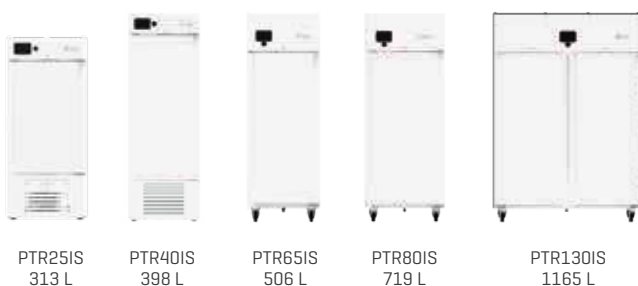This series offers high performance refrigerators for the storage of pharmaceutical preparations to meet the most demanding applications. In order to achieve that purpose, these models have incorporated **INVERTER** technology with high energy efficiency and low energy consumption and lower noise level. These models are designed according to **DIN 58435** standard. Capacity ranges from 313 to 1202 litres, and all models with Digital Plus. controller. Solid or glass doors are available in the whole range.

Pág. 30 - 33

CARACTERÍSTICAS GENERALES / GENERAL FEATURES

CONTROLADOR DIGITAL (modelos gama Pharma CARE) o CONTROLADOR DIGITAL PLUS (modelos gama Pharma CARE PLUS)

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DIGITAL CONTROLLER (Pharma CARE range models) or DIGITAL PLUS CONTROLLER (Pharma CARE and Pharma CARE PLUS range models)

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LOCK / DOOR SENSOR

All our equipment has locks and open door sensors, to guarantee the security of the stored products by controlling opening time and fans functioning.

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LED LIGHTING

The use of LED lighting reduces the energy consumption and improves the homogeneity of the temperature inside the cabinet.

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DOOR STYLE

It is possible to choose solid or glass doors in our whole range. All our doors are hinged and reversible.

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EASY REMOVABLE DOOR GASKET

It is ideal to maintain optimal hygiene and facilitate maintenance work.

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ROUNDED CORNERS

The interiors have rounded corners, which facilitate their cleaning.

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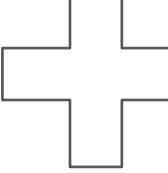
FLEXIBLE STORAGE SOLUTIONS

Our new shelf / drawer system allows you to adjust their position, vary their number, change shelves for drawers and vice versa or complement them with divisors. This provides the fit of the equipment to your needs and storage convenience of medications or vaccines of different sizes.

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BATTERY

Allows the maintenance of temperature and alarms register in case of a power failure.

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INVERTER

INVERTER technology (Pharma CARE PLUS range models) with high energy efficiency and low energy consumption.

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FREEZING PROTECTION ELECTRONIC CONTROLLER

Safety thermostat prevents the temperature dropping below +2°C.

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EVAPORATION PAN

Our equipment has a built-in tray to evaporate automatically defrosted water.

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ADVANCED DEFROST SYSTEM

Built-in to ensure uniform cooling over time and to facilitate maintenance. In addition, the equipment analyses its operation and only defrosts when it is necessary, to improve the stability of the temperature.

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CASTERS / ADJUSTABLE LEGS / HEIGHT REGULATORS

In our whole range, there is a possibility to incorporate casters to facilitate displacement and cleaning.

ARMARIO DE REFRIGERACIÓN PHARMA CARE / PHARMA CARE REFRIGERATOR



DIN 13277

GAMA SERIES	Pharma CARE	Pharma CARE	Pharma CARE	Pharma CARE	Pharma CARE	Pharma CARE
MODELO MODEL	PER07S	PER07G	PER16S	PER16G	PER24S	PER24G
Rango T° Temperature range [°C]	+2 / +8	+2 / +8	+2 / +8	+2 / +8	+2 / +8	+2 / +8
Capacidad [litros] Capacity [litres]	58	60	141	141	236	238
Medidas Exterior [mm] L x P x A Exterior dimensions [mm] W x D x H	445 x 526 x 818	445 x 508 x 818	600 x 688 x 824	600 x 680 x 824	600 x 688 x 1244	600 x 680 x 1244
Medidas Interior [mm] L x P x A Interior dimensions [mm] W x D x H	358 x 312 x 523	358 x 323 x 523	505 x 451 x 617	505 x 455 x 617	505 x 451 x 1037	505 x 455 x 1037
Claro de puerta [mm] L x A Door opening [mm] W x H	358 x 516	358 x 516	505 x 552	505 x 552	505 x 972	505 x 972
Espesor Aislamiento [mm] Insulation thickness [mm]	40	40	45	45	45	45
Acabado Exterior Exterior finishing	Acero galvanizado plastificado blanco White plastic coated galvanised steel					
Acabado Interior Interior finishing	Acero galvanizado plastificado blanco White plastic coated galvanised steel					
Nº Estantes Shelves Number	2	2	3	3	4	4
Dimensiones Estante [mm] L x P Shelf Size [mm] W x D	315 x 248	315 x 248	458 x 398	458 x 398	458 x 398	458 x 398
Material Estantes Shelves Material	Acero galvanizado con pintura Epoxi Epoxy coated galvanised steel					
Intercambiabilidad de Estantes/Cajones Interchangeable Shelves/Drawers	No	No	Si Yes	Si Yes	Si Yes	Si Yes
Tipo de Puerta Door Type	Ciega Solid	Cristal Glass	Ciega Solid	Cristal Glass	Ciega Solid	Cristal Glass
Cerradura Lock	Si Yes					
Salida USB USB Port	No					
Datalogger Datalogger	Opcional Optional					
Batería Battery	Si Yes					
Sonda simulación producto Product temperature simulation probe	Opcional Optional					
Sonda de temperatura ambiente Interior room temperature probe	Opcional Optional					
Termostato anticongelación Freezing protection controller	Si Yes					
Tipo de Controlador Controller type	Digital					
Tipo de Refrigerante Refrigerant type	R600a	R600a	R600a	R600a	R600a	R600a
Potencia Eléctrica [W] Electric power [W]	113	113	115	115	115	115
Tensión/Frecuencia [V / Hz] Voltage/Frequency [V / Hz]	230 / 50					
Peso [kg] Weight [kg]	45	49	62	65	92	96


* Pruebas realizadas con una temperatura ambiente de +25°C * Measured at +25°C ambient temperature

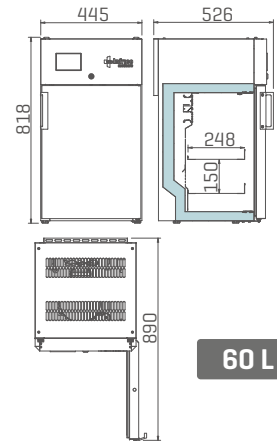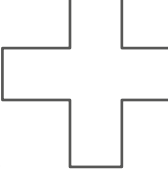Nos reservamos el derecho a modificar las especificaciones técnicas y constructivas sin previo aviso.
We reserve the right to modify technical and construction specifications without prior notice.


60 L

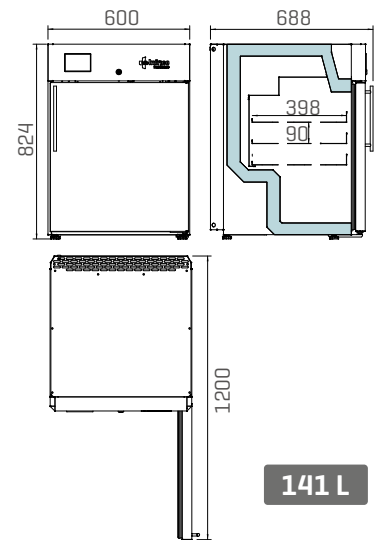PER07

141 L

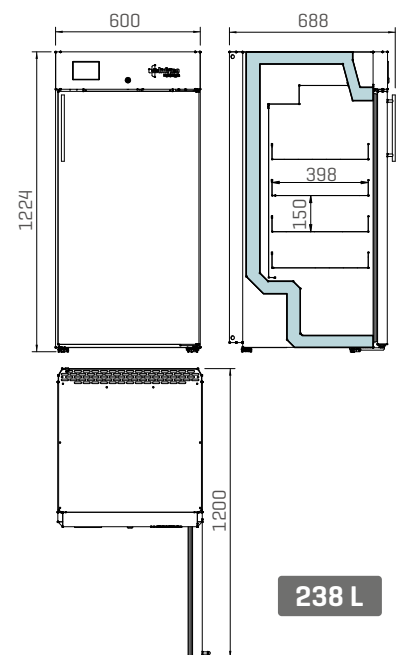PER16

238 L

PER24

ARMARIO DE REFRIGERACIÓN PHARMA CARE / PHARMA CARE REFRIGERATOR



DIN 13277

GAMA SERIES	Pharma CARE	Pharma CARE	Pharma CARE	Pharma CARE
MODELO MODEL	PER28S	PER28G	PER37S	PER37G
Rango T ^a Temperature range [°C]	+2 / +8	+2 / +8	+2 / +8	+2 / +8
Capacidad [litros] Capacity [litres]	327	330	373	376
Medidas Exterior [mm] L x P x A Exterior dimensions [mm] W x D x H	600 x 688 x 1644	600 x 680 x 1644	600 x 688 x 1844	600 x 680 x 1844
Medidas Interior [mm] L x P x A Interior dimensions [mm] W x D x H	505 x 451 x 1437	505 x 455 x 1437	505 x 451 x 1637	505 x 455 x 1637
Claro de puerta [mm] L x A Door opening [mm] W x H	505 x 1372	505 x 1372	505 x 1572	505 x 1572
Espesor Aislamiento [mm] Insulation thickness [mm]	45	45	45	45
Acabado Exterior Exterior finishing	Acero galvanizado plastificado blanco White plastic coated galvanised steel			
Acabado Interior Interior finishing	Acero galvanizado plastificado blanco White plastic coated galvanised steel			
Nº Estantes Shelves Number	5	5	6	6
Dimensiones Estante [mm] L x P Shelf Size [mm] W x D	458 x 398	458 x 398	458 x 398	458 x 398
Material Estantes Shelves Material	Acero galvanizado con pintura Epoxi Epoxy coated galvanised steel			
Intercambiabilidad de Estantes/Cajones Interchangeable Shelves/Drawers	Si Yes	Si Yes	Si Yes	Si Yes
Tipo de Puerta Door Type	Ciega Solid	Cristal Glass	Ciega Solid	Cristal Glass
Cerradura Lock	Si Yes			
Salida USB USB Port	No			
Datalogger Datalogger	Opcional Optional			
Batería Battery	Si Yes			
Sonda simulación producto Product temperature simulation probe	Opcional Optional			
Sonda de temperatura ambiente Interior room temperature probe	Opcional Optional			
Termostato anticongelación Freezing protection controller	Si Yes			
Tipo de Controlador Controller type	Digital			
Tipo de Refrigerante Refrigerant type	R290	R290	R290	R290
Potencia Eléctrica [W] Electric power [W]	210	210	210	210
Tensión/Frecuencia [V / Hz] Voltage/Frequency [V / Hz]	230 / 50			
Peso [kg] Weight [kg]	104	112	127	138


* Pruebas realizadas con una temperatura ambiente de +25°C * Measured at +25°C ambient temperature

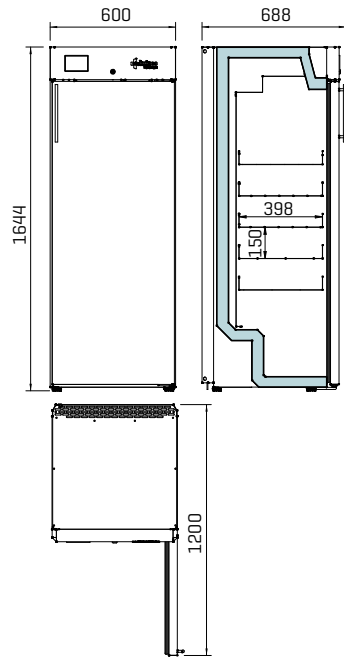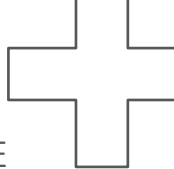Nos reservamos el derecho a modificar las especificaciones técnicas y constructivas sin previo aviso.
We reserve the right to modify technical and construction specifications without prior notice.

330L

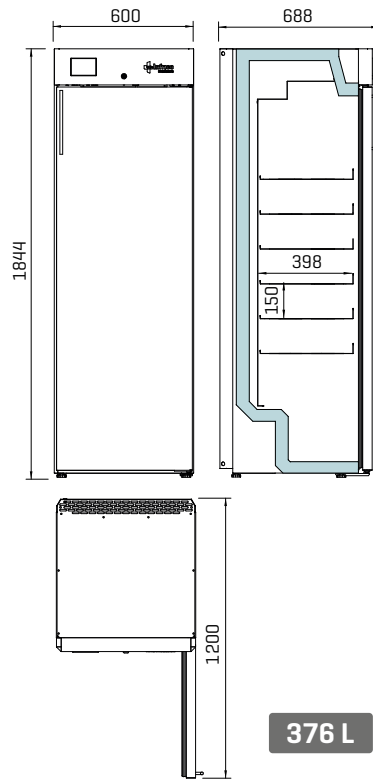PER28

376 L

PER37

ARMARIO DE REFRIGERACIÓN PHARMA CARE PLUS / PHARMA CARE PLUS REFRIGERATOR

DIN 13277


inverter


GAMA SERIES	Pharma CARE PLUS	Pharma CARE PLUS	Pharma CARE PLUS	Pharma CARE PLUS	Pharma CARE PLUS	Pharma CARE PLUS
MODELO MODEL	PTR25IS	PTR25IG	PTR40IS	PTR40IG	PTR65IS	PTR65IG
Rango T° Temperature range [°C]	+2 / +8	+2 / +8	+2 / +8	+2 / +8	+2 / +8	+2 / +8
Capacidad [litros] Capacity [litres]	251	261	398	413	506	521
Medidas Exterior [mm] L x P x A Exterior dimensions [mm] W x D x H	670 x 697 x 1538	670 x 697 x 1538	670 x 683 x 1963	670 x 683 x 1963	687 x 870 x 1955	687 x 870 x 1955
Medidas Interior [mm] L x P x A Interior dimensions [mm] W x D x H	568 x 484 x 913	568 x 504 x 913	568 x 484 x 1338	568 x 504 x 1338	534 x 633 x 1207	534 x 654 x 1207
Claro de puerta [mm] L x A Door opening [mm] W x H	568 x 855	568 x 855	568 x 1280	568 x 1280	534 x 1350	534 x 1350
Espesor Aislamiento [mm] Insulation thickness [mm]	50	50	50	50	75	75
Acabado Exterior Exterior finishing	Acero galvanizado plastificado blanco White plastic coated galvanised steel					
Acabado Interior Interior finishing	Acero galvanizado plastificado blanco White plastic coated galvanised steel			Acero Inoxidable Stainless steel		
Nº Cajones Drawers Number	4	4	7	7	7	7
Dimensiones Cajón [mm] L x P Drawer Size [mm] W x D	500 X 443	500 X 443	500 x 443	500 x 443	452 x 585	452 x 585
Material Cajones Drawers Material	Acero galvanizado con pintura Epoxi Epoxy coated galvanised steel			Acero Inoxidable Stainless steel		
Intercambiabilidad de Estantes/Cajones Interchangeable Shelves/Drawers	Si Yes	Si Yes	Si Yes	Si Yes	Si Yes	Si Yes
Tipo de Puerta Door Type	Ciega Solid	Cristal Glass	Ciega Solid	Cristal Glass	Ciega Solid	Cristal Glass
Cerradura Lock	Si Yes					
Salida USB USB Port	Si Yes					
Datalogger Datalogger	Si Yes					
Batería Battery	Si Yes					
Sonda simulación producto Product temperature simulation probe	Opcional Optional					
Sonda de temperatura ambiente Interior room temperature probe	Si Yes					
Termostato anticongelación Freezing protection controller	Si Yes					
Tipo de Controlador Controller type	Digital Plus					
Tipo de Refrigerante Refrigerant type	R600a	R600a	R600a	R600a	R290	R290
Potencia Eléctrica [W] Electric power [W]	133	138	133	152	426	445
Tensión/Frecuencia [V / Hz] Voltage/Frequency [V / Hz]	230 / 50					
Peso [kg] Weight [kg]	127	132	139	153	153	160

* Pruebas realizadas con una temperatura ambiente de +25°C
* Measured at +25°C ambient temperature

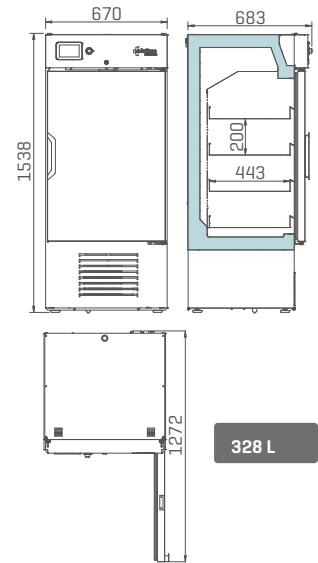
Nos reservamos el derecho a modificar las especificaciones técnicas y constructivas sin previo aviso.
We reserve the right to modify technical and construction specifications without prior notice.

PTR25

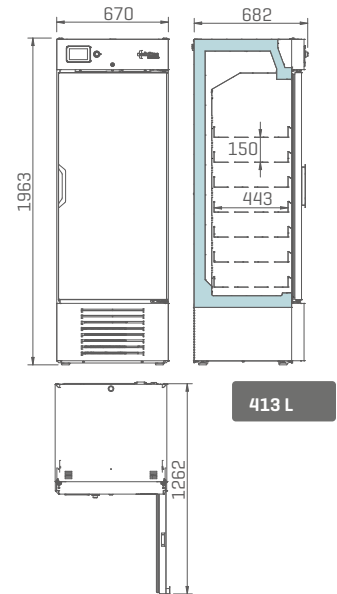PTR40


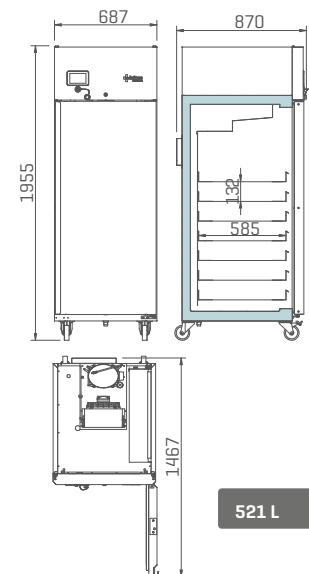
PTR65

DIN 13277


GAMA SERIES	Pharma CARE PLUS	Pharma CARE PLUS	Pharma CARE PLUS	Pharma CARE PLUS
MODELO MODEL	PTR80IS	PTR80IG	PTR130IS	PTR130IG
Rango T° Temperature range [°C]	+2 / +8	+2 / +8	+2 / +8	+2 / +8
Capacidad [litros] Capacity [litres]	719	737	1166	1202
Medidas Exterior [mm] L x P x A Exterior dimensions [mm] W x D x H	797 x 1000 x 1955	797 x 1000 x 1955	1385 x 834 x 1955	1385 x 834 x 1955
Medidas Interior [mm] L x P x A Interior dimensions [mm] W x D x H	645 x 803 x 1388	645 x 823 x 1388	1234 x 681 x 1388	1234 x 702 x 1388
Claro de puerta [mm] L x A Door opening [mm] W x H	645 x 1350	645 x 1350	534 x 1350	534 x 1350
Espesor Aislamiento [mm] Insulation thickness [mm]	75			
Acabado Exterior Exterior finishing	Acero galvanizado plastificado blanco White plastic coated galvanised steel			
Acabado Interior Interior finishing	Acero Inoxidable Stainless steel			
Nº Cajones Drawers Number	7	7	14	14
Dimensiones Cajón [mm] L x P Drawer Size [mm] W x D	564 x 750	564 x 750	452 x 585	452 x 585
Material Cajones Drawers Material	Acero Inoxidable Stainless steel			
Intercambiabilidad de Estantes/Cajones Interchangeable Shelves/Drawers	Si Yes			
Tipo de Puerta Door Type	Ciega Solid	Cristal Glass	Ciega Solid	Cristal Glass
Cerradura Lock	Si Yes			
Salida USB USB Port	Si Yes			
Datalogger Datalogger	Si Yes			
Batería Battery	Si Yes			
Sonda simulación producto Product temperature simulation probe	Opcional Optional			
Sonda de temperatura ambiente Interior room temperature probe	Si Yes			
Termostato anticongelación Freezing protection controller	Si Yes			
Tipo de Controlador Controller type	Digital Plus			
Tipo de Refrigerante Refrigerant type	R290			
Potencia Eléctrica [W] Electric power [W]	434	442	718	750
Tensión/Frecuencia [V / Hz] Voltage/Frequency [V / Hz]	230 / 50			
Peso [kg] Weight [kg]	172	178	208	225

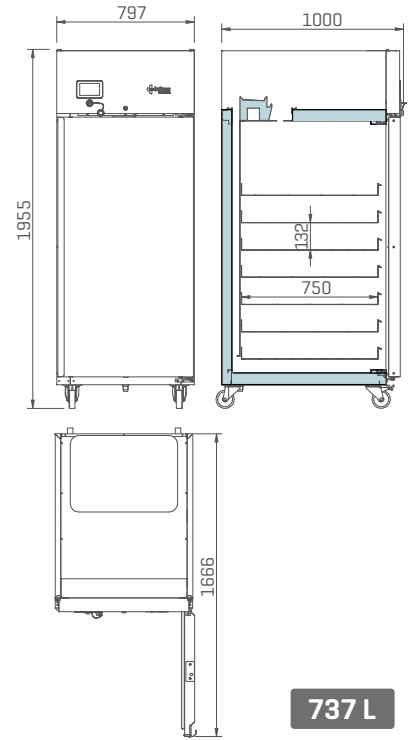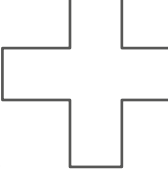* Pruebas realizadas con una temperatura ambiente de +25°C * Measured at +25°C ambient temperature

737 L

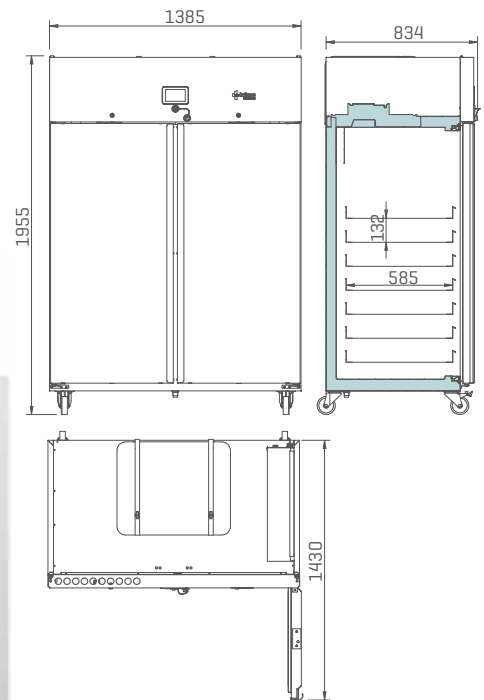PTR80I

1202 L

PTR130I

Distribuye:

Akralab, S.L.U.

Pol. Ind. Las Atalayas
Av. de la Antigua Peseta, 77
03114 Alicante (Spain)

atencion.clientes@akralab.es

+34 965 116 521

www.akralab.es